

SD 06 Insert cup cpl. For ultrasonic bath



Glass, 600 ml, for insertion into perforated lid/insertion strip with ring GR 06 and lid DD 06. Insertion beakers are used when applying chemically aggressive liquids (e.g. TICKOPUR J 80 U), removing chemically aggressive or abrasive contaminants or sonicating sample liquids (indirect sonication). Because nothing may be placed on the bottom of the ultrasonic bath, BANDELIN offers suitable perforated lids/insert strips to hold the insert beakers. Further advantage: Several insert cups can be sonicated with different liquids at the same time (e.g. removal of dental cements with STAMMOPUR Z in one and cleaning with simultaneous chemical disinfection of dental instruments with STAMMOPUR DR 8 in a second insert cup). Please note that in order to transfer the ultrasound to the insert cup, the surface tension of the water must be reduced and therefore a cleaning solution (e.g. TICKOPUR R 33 3%) must be used as coupling liquid / contact liquid.

- for hanging in perforated cover/insert strip
- suitable for various SONOREX ultrasonic baths
- D = 84 mm, H = 125 mm
- capacity = 600 ml

- Glass, complete with retaining ring and lid.
- for indirect cleaning of small parts

Item no.	WL17905
Model	SD 06
Manufacturer	BANDELIN
Manufacturer article no.	330
GPSR manufacturer data	BANDELIN electronic GmbH & Co. KG Heinrichstrasse 3-4 DE-12207 Berlin www.bandelin.com
Length	84 mm
Width	90 mm
Height	125 mm
Sales unit	1 piece
Content unit	1 piece

