

PD 06 Plastic cup cpl. for ultrasonic bath

Plastic, 600ml, to hang in perforated lid DE. Use when using a cleaner or solvent that is aggressive for stainless steel (e.g. TICKOPUR J 80 U) is necessary to protect the cleaning device. Because the insert vessels must not be placed on the bottom of the tank, BANDELIN offers suitable perforated lids to hold the insert vessels. Another advantage of using several insert vessels in one ultrasonic unit is the possibility of using different cleaning and disinfecting preparations at the same time without changing the bath. (e.g. removal of dental cements with STAMMOPUR Z and disinfection and cleaning of instruments with STAMMOPUR DR 8 in the dental practice). Please note that in order to transfer the ultrasound to the insert vessel, the surface tension of the water must be reduced and therefore a cleaning solution (e.g. TICKOPUR R 33 3%) must be used as coupling liquid / contact liquid.

- for hanging in perforated cover/insert strip
- suitable for various SONOREX ultrasonic baths
- D = 86 mm, H = 125 mm
- capacity = 600 ml
- plastic, complete with retaining ring and lid
- for indirect cleaning of small parts

item number	WL18296
model	PD 06
manufacturer	BANDELIN
manufacturer item number	299
GPSR manufacturer information	BANDELIN electronic GmbH & Co. KG Heinrichstrasse 3-4 DE-12207 Berlin www.bandelin.com
length	86 mm
width	90 mm
height	125 mm
order unit	1 piece
content unit	1 piece